


## 【USB Driver Installation Procedure】 for Windows10

**Two types of USB drivers** are required to be installed before connecting Lumitester PD-30 and PC.  
Please follow the procedure below.

1 Connect Lumitester PD-30 to PC with the USB cable provided.


### 2 Install the first USB driver.

Click "Start" button and open the "Connected devices".  
Then Click "Device manager".

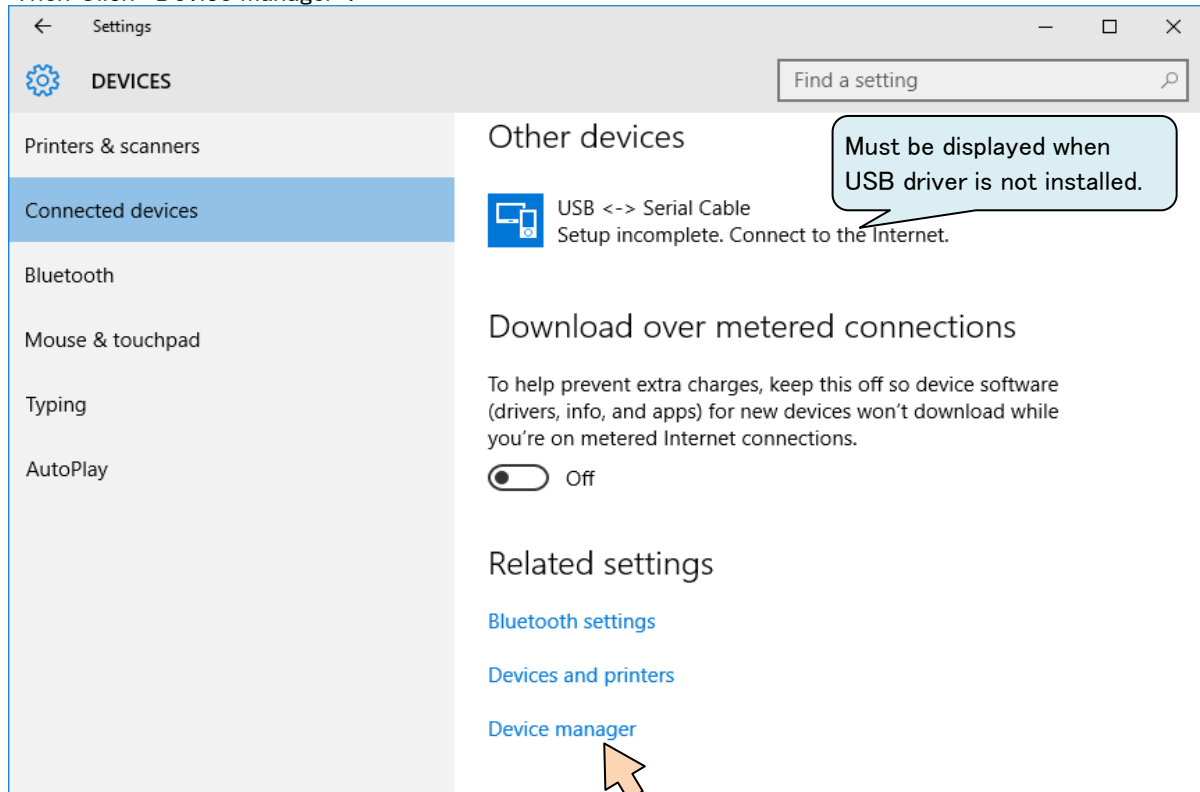


3 The device with the "⚠️" will be displayed on Device Manager.


Place the cursor on the name of device with "⚠️" and right-click, and select "Update Driver Software".

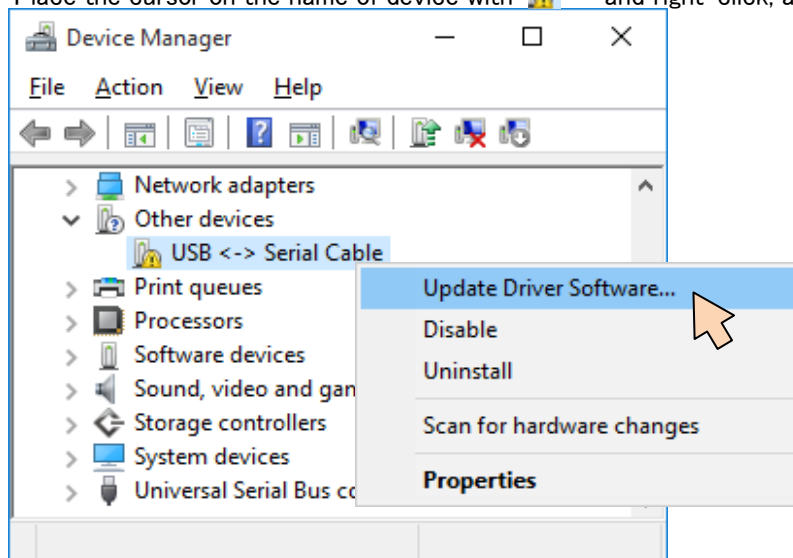


- 4 "Update Driver Software" is displayed on the screen.  
Click "Browse my computer for driver software".

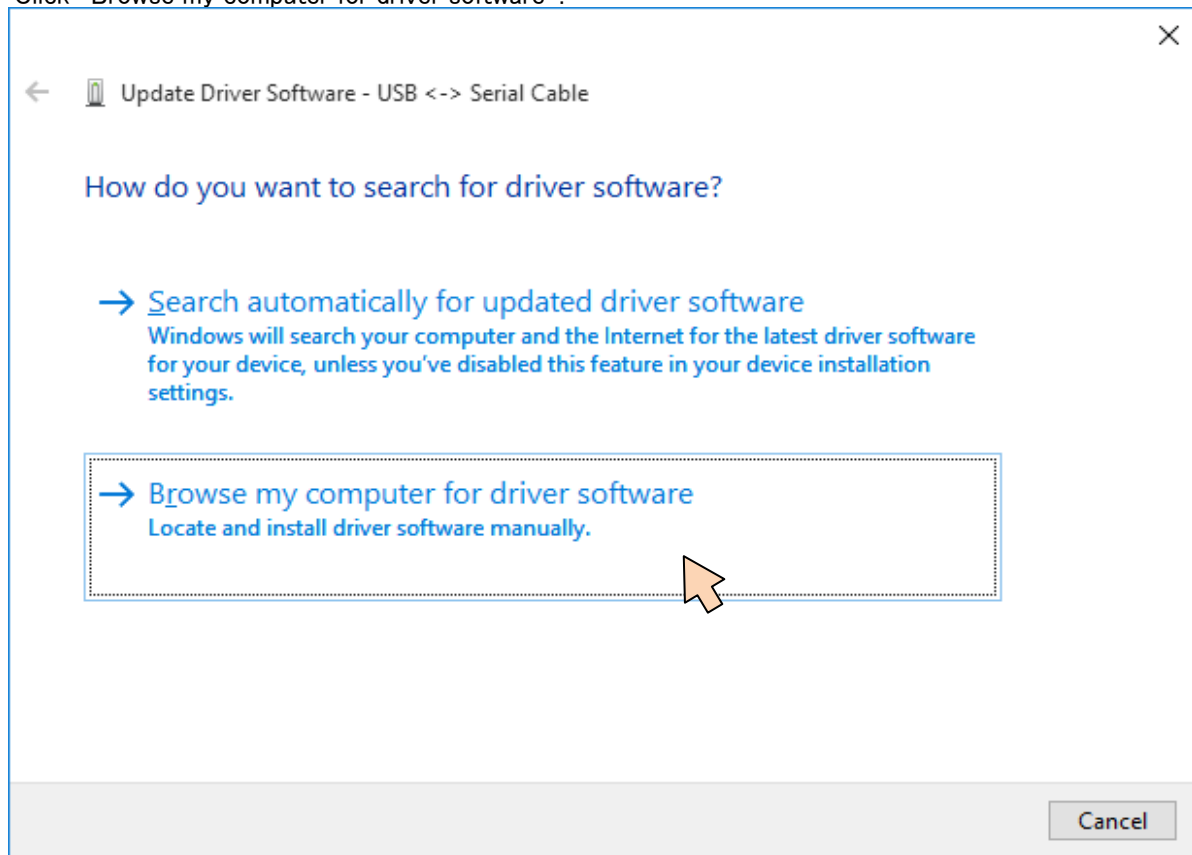

- 5 Click "Browse", find and select the folder of "USB Drivers" downloaded, and then click "Next".

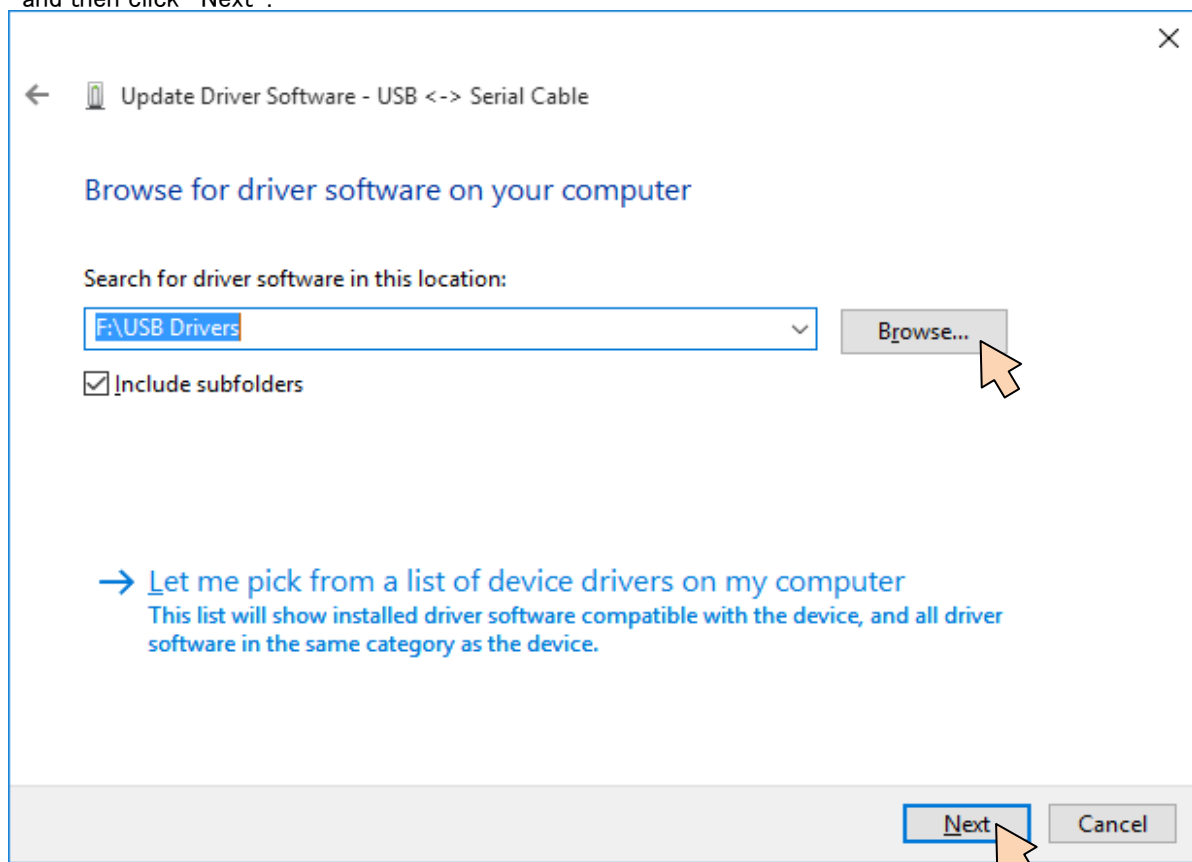


6 Installation of the driver software will begin.

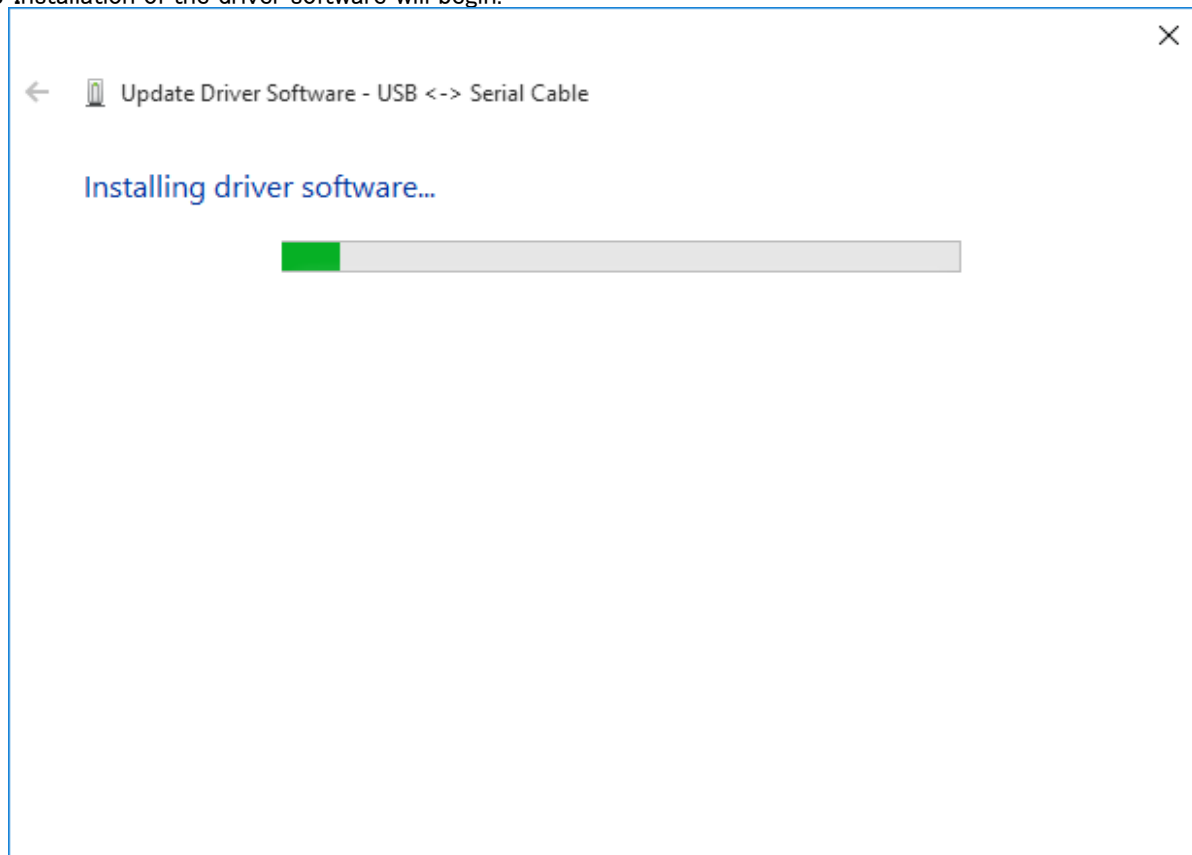


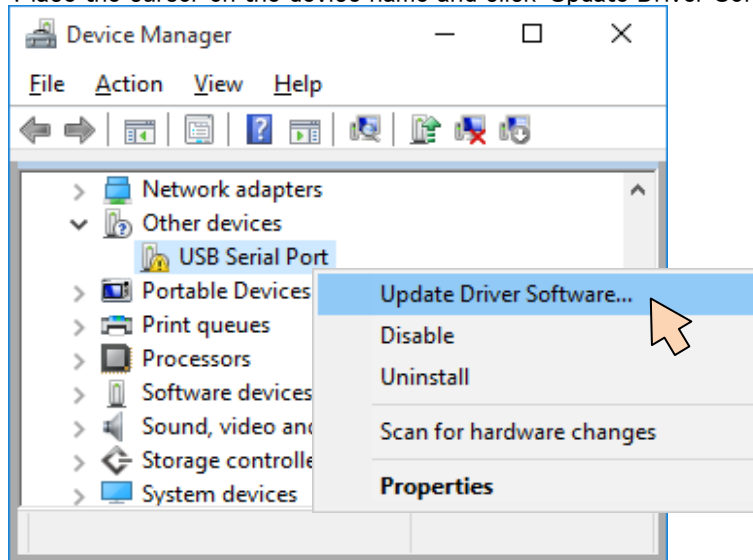
7 The **first USB driver** is now installed.

Click "Close" to close the window. And begin installation for the **second USB driver**.

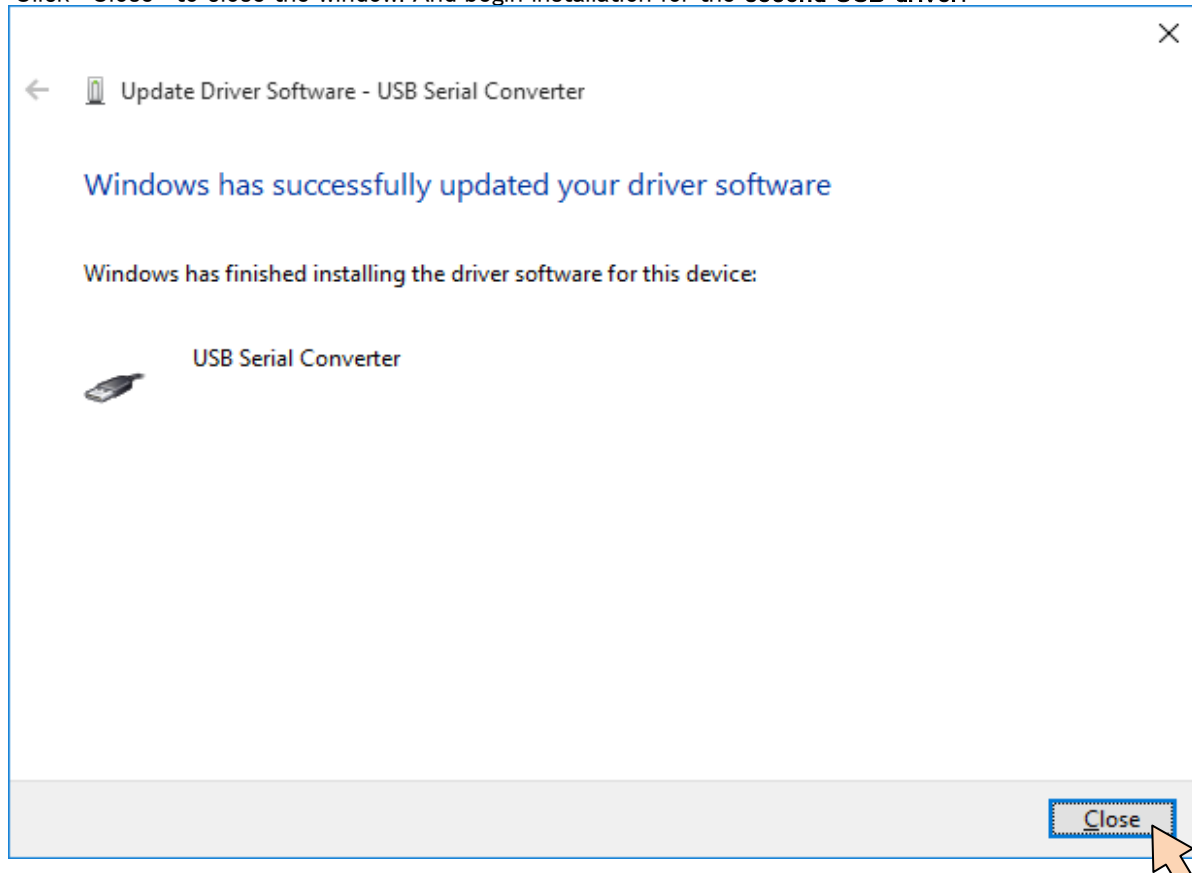

8 For the Installation of **second USB driver**.

Open the Device Manager in the procedure. The device name with the  " is displayed on Device Manager. Place the cursor on the device name and click "Update Driver Software".



9 Follow the same way as in the procedure 4 to 7 of first USB driver for installation.

When the installation is completed the following screen appears.

Click "Close" to close the window.

